Emergency Eyewash Station Weekly Inspection Checklist

How to Inspect & Test:

- Is the area around the station clear of obstructions?
- Is the eyewash clean and free of dirt/debris?
- Flush eyewash for 30 seconds weekly
- Do the covers/caps open without assistance?
- Is the plumbing free from leakage?
- Are the jets working properly (both spigots are delivering evenly, and with sufficient but not too great of a flow)?
- Is the water clean and clear?

How to Obtain Assistance:

Questions, maintenance, training EHS: 302-831-8475 | dehsafety@udel.edu

Maintenance

requests may also be called in directly **Facilities Services**: 302-831-1141

Emergencies:

Hold eyes open and flush for at least 15 minutes. CALL 911 and seek medical assistance

	as the water death and clour.						
Date	Initials	Comment		Date	Initials	Comment	